

EPA 3051 (for oil)

Milestone Application Note for Acid Digestion SK-Officials-002



Summary

This method represents an Official method for the microwave acid digestion of Oil samples with the Milestone ETHOS UP and ETHOS EASY systems.

Instrumentation

Milestone ETHOS UP or ETHOS EASY microwave units, SK-15 high-pressure rotor, T1 temperature control. Options T2 temperature control and P1 or P2 pressure control.



Sample weight and reagents

Sample weight	Up to 250 mg
Reagents	10 mL HNO ₃

Microwave program

Step	Time	T1	Power
1	00:10:00	180°C	1800 W*
2	00:10:00	180°C	1800 W*

Subject to change without notice.
For additional information please contact
application@milestonesrl.com

General precautions

*Use only up to 500 W for runs with 3 or less vessels simultaneously.

Always use hand, eye and body protection when operating with the microwave system.

