

ASTM D1506 (Method B)

Milestone Application Note for Ashing – High Throughput - Officials



Summary

This procedure covers the determination of the ash content of carbon black samples according to ASTM D1506 (B) method.

Instrumentation

Milestone PYRO microwave system


HighThroughput muffle furnace

Crucibles (ceramic, quartz, quartz fibre)

Analytical balance ($\pm 0,0001$ g)

Disseccator

Spatula

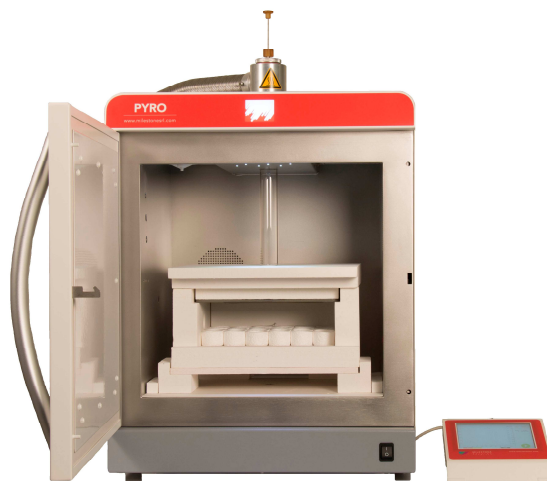

Sample amount

| | |
|---------------|-----|
| Sample weight | 2 g |
|---------------|-----|

Microwave program

| Step | Time | T1 | Power | Vent. |
|------|----------|-------|--------|-------|
| 1 | 00:25:00 | 550°C | 1800 W | * |
| 2 | 01:00:00 | 550°C | 1800 W | * |

Procedure

1. Place a clean crucible on the balance plate, tare it and weigh the sample.
2. Place the crucible into the preheated furnace.
3. Remove the crucible from the furnace and cool it in dissector.
4. Place the crucible on the balance plate and take the weigh.
5. Calculate the ash residue.

Notes

The crucible before use must be at constant weight.

All modifications to the present method must be agreed with Milestone Application Specialist.

General precautions

Always use hand, eye and body protection when operating with the microwave system.

Subject to change without notice.
For additional information please contact
application@milestonesrl.com

